

# In-Fusion® SMARTer® Directional Cloning & Screening Components

Catalog No.

Lot Number

639632 (Not sold separately)

Specified on product label.

#### **Product Information**

This package contains the components needed for cloning In-Fusion SMARTer Directional cDNA Libraries.

#### **Package Contents**

- 60 μl pSMART2IFD Linearized Vector (150 ng/μl)
- 100 μl Forward Screening Primer (10 μM)
- 100 μl Reverse Screening Primer (10 μM)

#### **Storage Conditions**

Store all components at -20°C.

#### **Shelf Life**

1 year from date of receipt under proper storage conditions.

### **Shipping Conditions**

Dry ice  $(-70^{\circ}\text{C})$ 

## pSMART2IFD Linearized Vector

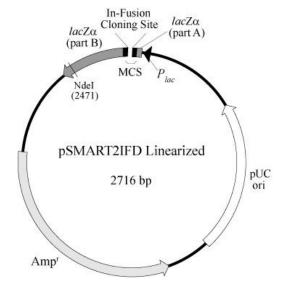


Figure 1. pSMART2IFD Linearized Vector Map

Clontech Laboratories, Inc.

A Takara Bio Company

1290 Terra Bella Avenue, Mountain View, CA 94043, USA


U.S. Technical Support: tech@clontech.com

United States/Canada Asia Pacific 800.662.2566 (012513)

+1.650.919.7300

Europe +33.(0)1.3904.6880

Japan +81.(0)77.543.6116 In-Fusion® SMARTer® Directional Cloning & Screening Components (Not sold separately)

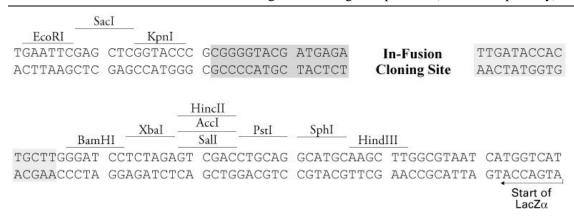

**Figure 2.** pSMART2IFD Linearized Vector In-Fusion Cloning Site. Each shaded region contains 15 bp of sequence that is complementary to one of the ends of the cDNA generated with the In-Fusion SMARTer PCR cDNA Synthesis Kit. The ends are different to allow for directional cDNA cloning.

## **Description**

The pSMART2IFD Linearized Vector is a pUC19-based, high copy number, *E. coli* cloning vector. This linearized vector was generated by PCR, and contains the blunt ends shown flanking the In-Fusion Cloning Site in the sequence above. Each end of the vector shares 15 bp of complementarity (shaded above) with a different end of the cDNA generated using the In-Fusion SMARTer PCR cDNA Synthesis Kit (Cat. No. 634934). The ends are different to allow for directional cDNA cloning.

The In-Fusion Cloning Site is located within the  $lacZ\alpha$  gene (encoding the N-terminal fragment of  $\beta$ -galactosidase), the expression of which is driven by the lac promoter ( $P_{lac}$ ). The location of the cloning site within  $lacZ\alpha$  allows for blue/white selection (i.e.,  $\alpha$ -complementation) after In-Fusion cloning. The vector also contains a pUC origin of replication that allows propagation in E. coli and temperature-dependent regulation of vector copy number, and an ampicillin resistance gene (Amp<sup>r</sup>) for selection in E. coli.

#### **Location of Features**

- $lacZ\alpha$  (N-terminal fragment of  $\beta$ -galactosidase; part A): 1–69 (complementary)
- $P_{lac}$  (lac promoter): 114–143 (complementary)
- pUC origin of replication: 467–1055 (complementary)
- Amp<sup>r</sup> (ampicillin resistance gene; β-lactamase): 1226–2086 (complementary)
- lacZα (N-terminal fragment of β-galactosidase; part B): 2432–2716 (complementary)

#### **Additional Information**

The pSMART2IFD Linearized Vector is provided as part of the In-Fusion SMARTer Directional cDNA Library Construction Kit (Cat. No. 634933), and is designed for effortless library construction with In-Fusion cloning technology.

#### Propagation in *E. coli*

- Recommended host strain: DH5α, XL1-Blue, and other general purpose strains.
- Selectable marker: plasmid confers resistance to ampicillin (50 μg/ml) in E. coli hosts.
- E. coli replication origin: pUC
- Copy number: high
- Cell transformation efficiency: > 4 x 10<sup>9</sup> cfu/μg

(012513) Page 2 of 3

In-Fusion® SMARTer® Directional Cloning & Screening Components (Not sold separately)

#### References

1. Yanisch-Perron, C., et al. (1985) Gene 33(1):103-119.

## **Quality Control Data**

Double-stranded SMARTer cDNA was synthesized from Control Mouse Liver Total RNA as described in the In-Fusion SMARTer Directional cDNA Library Construction Kit User Manual (PT5147-1). Three different volumes of SMARTer cDNA were each combined with 2  $\mu$ l (150 ng/ $\mu$ l) of pSMART2IFD in separate In-Fusion cloning reactions, according to the user manual. Each cloning reaction was transformed into a culture of *E. coli*. The three test reactions yielded at least  $1x10^6$  independent colonies, of which  $\geq 75\%$  were white. Out of 15 randomly selected white colonies screened by PCR, 10 contained inserts  $\geq 700$  bp in length.

#### **Plasmid Identity & Purity**

Vector identity was confirmed by sequencing.

•  $A_{260}/A_{280}$ : 1.8–2.0

(012513) Page 3 of 3

## **Notice to Purchaser**



# In-Fusion® SMARTer® Directional Cloning & Screening Components

CATALOG NO.

639632

#### NOTICE TO PURCHASER:

Our products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, in vitro diagnostic purposes, therapeutics, or in humans. Our products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without prior written approval of Clontech Laboratories, Inc.

Your use of this product is also subject to compliance with the licensing requirements listed below and described on the product's web page at <a href="http://www.clontech.com">http://www.clontech.com</a>. It is your responsibility to review, understand and adhere to any restrictions imposed by these statements.

#### TRADEMARKS:

Clontech®, the Clontech logo, In-Fusion®, and SMARTer® are trademarks of Clontech Laboratories, Inc.

All other trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech is a Takara Bio Company. ©2016 Clontech Laboratories, Inc. This document has been reviewed and approved by the Clontech Quality Assurance Department.

Clontech Laboratories, Inc.

A Takara Bio Company 1290 Terra Bella Avenue, Mountain View, CA 94043, USA U.S. Technical Support: tech@clontech.com

3/30/2016

United States/Canada Asia Pacific Europe Japan

800.662.2566 +1.650.919.7300 +33.(0)1.3904.6880 +81.(0)77.543.6116