

I. List of Components

Store all components at -20°C .

- MMLV Reverse Transcriptase, GPR
- 5X Reverse Transcription Buffer
- 100 mM DTT

II. Additional Materials Required

- RNase-free H_2O
- dNTP Mix (10 mM each); we recommend Advantage® UltraPure PCR Deoxynucleotide Mix (Cat. No. 639125)
- Oligo (dT)₁₂₋₁₈, random hexamer or gene-specific primers may be used.
- [Optional] 60 mM EDTA

III. For Routine First-Strand cDNA Synthesis Reactions

This 20- μl reaction is suitable for synthesizing first-strand cDNA from 5 ng–5 μg of total RNA or 10 ng–1 μg of poly A⁺ RNA.

1. Add 2.5 μl of 20 μM primer stock (final concentration 2.5 μM) to your RNA sample. Add RNase-free H_2O to a final volume of 11.5 μl .
2. Heat the mixture at 70°C for 3 min, and immediately cool on ice.
3. Centrifuge briefly, then add the following:
 - 4 μl 5X Reverse Transcription Buffer
 - 2 μl dNTP Mix
 - 2 μl 100 mM DTT
4. Mix the contents of the tube by gently pipetting up and down. Add 0.5 μl MMLV Reverse Transcriptase, GPR and mix again.
5. Incubate at 42°C for 60 min.

NOTE: Samples can be incubated for 50–90 min if necessary.

6. Terminate the reaction by heating at 70°C for 15 min, or by adding 4 μl of 60 mM EDTA.

IV. References

Ausubel, F. M., *et al.* (1995) In *Current Protocols in Molecular Biology* (Greene Publishing Associates, Inc. and John Wiley & Sons, Inc.), Supplement 29, Section 5.5.

Contact Us	
Customer Service/Ordering	Technical Support
Telephone: 800.662.2566 (toll-free)	Telephone: 800.662.2566 (toll-free)
Fax: 800.424.1350 (toll-free)	Fax: 650.424.1064
Web: www.clontech.com	Web: www.clontech.com
E-mail: orders@clontech.com	E-mail: tech@clontech.com

Notice to Purchaser

General Purpose Reagent. This product is intended For Laboratory Use. Outside of the United States, this product is intended for research use only unless otherwise stated. This product is not intended for a specific application or made for any clinical use. The performance characteristics of this product have not been fully established. It is the user's responsibility to validate the performance of the product, and any component thereof, for any particular use. Resale or transfer of this product, any component thereof, or any substance produced through use of this product, or any component thereof, to any third party is expressly forbidden. To obtain additional rights, please contact BD_OEM@clontech.com.

Your use of this product is also subject to compliance with the licensing requirements listed below and described on the product's web page at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by these statements.

Clontech®, the Clontech logo, and Advantage® are trademarks of Clontech Laboratories, Inc. All other trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech Laboratories, Inc. is a Takara Bio Company. ©2016 Clontech Laboratories, Inc.

This document has been reviewed and approved by the Clontech Quality Assurance Department.