

Titanium® *Taq* PCR Kit Protocol-at-a-Glance

(PT3304-2)

Please read the *User Manual* before using this Protocol-at-a-Glance. This abbreviated protocol is provided for your convenience, but is not intended for first-time users.

Primer Design

Primer design is the single largest variable in PCR applications and the single most important factor in determining the success or failure of PCR reactions. *Always check your primer design before constructing or ordering primers.*

Titanium *Taq* can be used in a wide variety of PCR applications, and the constraints on primer design will vary from one application to the next. In general, primers should have a T_m of $\sim 70^\circ\text{C}$ to achieve optimal results in a two-step cycling program with a 68°C combined annealing/extension step. Therefore, whenever possible, primers should be *at least* 22 nucleotides long (25–30-mers are preferred) and should have a GC content of 45–60%. Furthermore, the 3' ends of each primer should not be complementary to each other and should have a low GC content.

Setting up the Reaction

Combine the following in a PCR tube on ice:

Volume	Reagent
40 μl	PCR-Grade Water
5 μl	10X Titanium <i>Taq</i> PCR Buffer
1 μl	50X dNTP Mix (10 mM ea.)
1 μl	5' primer (10 μM)
1 μl	3' primer (10 μM)
1 μl	50X Titanium <i>Taq</i> DNA Polymerase
1 μl	DNA Template (100 ng/ μl)
50 μl	Total volume

Recommended Cycling Conditions

Use the following guidelines when setting up your initial experiments with Titanium *Taq*. These are general guidelines—the optimal cycling conditions may vary with different thermal cyclers and will depend on your particular primers, templates, and other experimental variables.

95°C for 1 min	} 1 cycle
95°C 30 sec	} 25–35 cycles
68°C 1–3 min	
68°C for 3 min	} 1 cycle



Clontech

United States/Canada
800.662.2566

Asia Pacific
+1.650.919.7300

Europe
+33.(0)1.3904.6880

Japan
+81.(0)77.543.6116

Clontech Laboratories, Inc.
A Takara Bio Company
1290 Terra Bella Ave.
Mountain View, CA 94043
Technical Support (US)
E-mail: tech@clontech.com
www.clontech.com

Recommendations for Electrophoresis

Expected insert size range	% agarose	Recommended DNA size markers
0.3–1.5 kb	1.5	$\phi\text{X174}/\text{Hae III}$
0.5–10 kb	1.2	1 kb DNA ladder
>5 kb	0.8	$\lambda/\text{Hind III}$

(PR163867; published June 2011)

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product's web page at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

Clontech, the Clontech logo, and Titanium are trademarks of Clontech Laboratories, Inc. All other marks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Clontech is a Takara Bio Company. ©2011 Clontech Laboratories, Inc.

This document has been reviewed and approved by the Clontech Quality Assurance Department.