



Restriction Map and Multiple Cloning Site (MCS) of pDsRed-Monomer-C1. Unique restriction sites are in bold. NOTE: The *Xba*I and *Bcl*I sites are methylated in the DNA provided by Clontech Laboratories, Inc. If you wish to digest the vector with these enzymes, you will need to transform the vector into a *dam*⁻ host and make fresh DNA.

Description

pDsRed-Monomer-C1 is a mammalian expression vector that encodes DsRed-Monomer (DsRed-M1), a monomeric mutant derived from the tetrameric *Discosoma* sp. red fluorescent protein DsRed (1). DsRed-Monomer contains a total of forty-five amino acid substitutions (listed on page 2). When DsRed-Monomer is expressed in mammalian cell cultures, red fluorescent cells can be detected by either fluorescence microscopy or flow cytometry 12–16 hr after transfection (DsRed-Monomer excitation and emission maxima = 557 nm and 592 nm, respectively). The DsRed-Monomer coding sequence is human codon-optimized for high expression levels in mammalian cells (2).

The multiple cloning site (MCS) in pDsRed-Monomer-C1 is positioned between the DsRed-Monomer coding sequence and the SV40 polyadenylation signal (SV40 poly A). Genes cloned into the MCS will be expressed as fusions to the C-terminus of DsRed-Monomer if they are in the same reading frame as DsRed-Monomer and there are no intervening stop codons. A Kozak consensus translation initiation site upstream of DsRed-Monomer increases the translation efficiency in eukaryotic cells (3). SV40 poly A signals downstream of the MCS direct proper processing of the 3' end of mRNA transcripts. The vector backbone also contains an SV40 origin for replication in mammalian cells expressing the SV40T-antigen, a pUC origin of replication for propagation in *E. coli*, and an f1 origin for single-stranded DNA production. A neomycin resistance cassette (Neo^r)—consisting of the SV40 early promoter, the neomycin/kanamycin resistance gene of Tn5, and polyadenylation signals from the Herpes simplex virus thymidine kinase (HSVTK) gene—allows stably transfected eukaryotic cells to be selected using G418. A bacterial promoter upstream of this cassette expresses kanamycin resistance in *E. coli*.



Clontech

United States/Canada
800.662.2566

Asia Pacific
+1.650.919.7300

Europe
+33.(0)1.3904.6880

Japan
+81.(0)77.543.6116

Clontech Laboratories, Inc.
A Takara Bio Company
1290 Terra Bella Ave.
Mountain View, CA 94043
Technical Support (US)
E-mail: tech@clontech.com
www.clontech.com

(PR0X3703; published October 2010)

Use

pDsRed-Monomer-C1 can be used to construct fusions to the C-terminus of DsRed-Monomer. If a fusion construct retains the fluorescent properties of the native DsRed-Monomer protein, its expression can be monitored by flow cytometry and its localization *in vivo* can be determined by fluorescence microscopy. The target gene must be cloned into pDsRed-Monomer-C1 so that it is in frame with the DsRed-Monomer coding sequences, with no intervening in-frame stop codons. The recombinant DsRed-Monomer vector can be transfected into mammalian cells using any standard transfection method. If required, stable transformants can be selected using G418 (4). pDsRed-Monomer-C1 can also be used as a cotransfection marker; the unmodified vector will express DsRed-Monomer.

We recommend using the DsRed-Monomer-C sequencing primer (see Sequencing primer location information below) to sequence genes cloned adjacent to the 3' end of the DsRed-Monomer coding region.

For Western blotting, the Living Colors® DsRed Polyclonal Antibody (Cat. No. 632496) can be used to recognize the DsRed-Monomer protein. However, for optimal results, it may be necessary to use a higher concentration of antibody than recommended on the DsRed Polyclonal Antibody Certificate of Analysis.

Location of features

- Human cytomegalovirus (CMV) immediate early promoter: 1–589
 - Enhancer region: 59–465; TATA box: 554–560
 - Transcription start point: 583
 - C→G mutation to remove *Sac* I site: 569
- Human codon-optimized DsRed-Monomer gene
 - Kozak consensus translation initiation site: 606–616
 - Start codon (ATG): 613–615; Stop codon: 1366–1368
 - Amino acid substitutions (DsRed→DsRed-Monomer)
 - GCC→GAC (Ala-2 to Asp) mutation: 616–618
 - TCC→AAC (Ser-3 to Asn) mutation: 619–621
 - TCC→ACC (Ser-4 to Thr) mutation: 622–624
 - AAG→GAG (Lys-5 to Glu) mutation: 625–627
 - AAC→GAC (Asn-6 to Asp) mutation: 628–630
 - CGC→CAG (Arg-13 to Gln) mutation: 649–651
 - ACC→TCC (Thr-21 to Ser) mutation: 673–675
 - GAG→TAC (Glu-26 to Tyr) mutation: 688–690
 - CGC→AAG (Arg-36 to Lys) mutation: 718–720
 - CAC→ACC (His-41 to Thr) mutation: 733–735
 - AAC→CAG (Asn-42 to Gln) mutation: 736–738
 - GTG→GCC (Val-44 to Ala) mutation: 742–744
 - AAG→CAG (Lys-47 to Gln) mutation: 751–753
 - GTG→GCC (Val-71 to Ala) mutation: 823–825
 - AAG→ATG (Lys-83 to Met) mutation: 859–861
 - AAG→ACC (Lys-92 to Thr) mutation: 886–888
 - GTG→TCC (Val-96 to Ser) mutation: 898–900
 - ACC→GAG (Thr-106 to Glu) mutation: 928–930
 - ACC→CAG (Thr-108 to Gln) mutation: 934–936
 - TCC→ACC (Ser-117 to Thr) mutation: 961–963
 - ATC→AAG (Ile-125 to Lys) mutation: 985–987
 - TCC→GCC (Ser-131 to Ala) mutation: 1003–1005
 - ATG→GCC (Met-141 to Ala) mutation: 1033–1035
 - GCC→CCC (Ala-145 to Pro) mutation: 1045–1047
 - CGC→AAG (Arg-149 to Lys) mutation: 1057–1059
 - CGC→CAG (Arg-153 to Gln) mutation: 1069–1071
 - CAC→TCC (His-162 to Ser) mutation: 1096–1098
 - AAG→CAC (Lys-163 to His) mutation: 1099–1101
 - CTG→ACC (Leu-174 to Thr) mutation: 1132–1134
 - GTG→TGC (Val-175 to Cys) mutation: 1135–1137
 - GAG→GAC (Glu-176 to Asp) mutation: 1138–1140
 - TCC→ACC (Ser-179 to Thr) mutation: 1147–1149
 - ATC→GTG (Ile-180 to Val) mutation: 1150–1152
 - ATG→AAG (Met-182 to Lys) mutation: 1156–1158
 - TAC→AAC (Tyr-192 to Asn) mutation: 1186–1188

TAC→CAC (Tyr-193 to His) mutation: 1189–1191
 TCC→AAC (Ser-203 to Asn) mutation: 1219–1221
 ATC→GTG (Ile-210 to Val) mutation: 1240–1242
 CGC→CAC (Arg-216 to His) mutation: 1258–1260
 ACC→GCC (Thr-217 to Ala) mutation: 1261–1263
 GGC→GCC (Gly-219 to Ala) mutation: 1267–1269
 CAC→TCC (His-222 to Ser) mutation: 1276–1278
 CTG→GGC (Leu-223 to Gly) mutation: 1279–1281
 TTC→TCC (Phe-224 to Ser) mutation: 1282–1284
 CTG→CAG (Leu-225 to Gln) mutation: 1285–1287

- Multiple Cloning Site: 1288–1368
- SV40 early mRNA polyadenylation signal
 Polyadenylation signals: 1507–1512 & 1536–1541; mRNA 3' ends: 1545 & 1657
- f1 single-strand DNA origin: 1604–2059 (Packages the noncoding strand of DsRed-Monomer)
- Bacterial promoter for expression of Kan^r gene
 –35 region: 2121–2126; –10 region: 2144–2149
 Transcription start point: 2156
- SV40 origin of replication: 2400–2535
- SV40 early promoter
 Enhancer (72-bp tandem repeats): 2233–2304 & 2305–2376
 21-bp repeats: 2380–3000, 2401–2421 & 2423–2440
 Early promoter element: 2456–2462
 Major transcription start points: 2452, 2490, 2496 & 2501
- Kanamycin/neomycin resistance gene
 Neomycin phosphotransferase coding sequences:
 Start codon (ATG): 2585–2587; stop codon: 3377–3379
 G→A mutation to remove *Pst* I site: 2767
 C→A (Arg to Ser) mutation to remove *Bss*H II site: 3113
- Herpes simplex virus (HSV) thymidine kinase (TK) polyadenylation signal
 Polyadenylation signals: 3615–3620 & 3628–3633
- pUC plasmid replication origin: 3964–4607

Sequencing primer location

- DsRed-Monomer-C sequencing primer (5'-AGCTGGACATCACCAACCACAACG-3'): 1205–1228
Note: The DsRed1-C Sequencing Primer (Cat. No. 632388) **cannot** be used as a sequencing primer for pDsRed-Monomer-C1.

Propagation in *E. coli*

- Suitable host strains: DH5 α , HB101, and other general purpose strains. Single-stranded DNA production requires a host containing an F plasmid such as JM109 or XL1-Blue.
- Selectable marker: plasmid confers resistance to kanamycin (50 μ g/ml) in *E. coli* hosts.
- *E. coli* replication origin: pUC
- Copy number: high
- Plasmid incompatibility group: pMB1/Col E1

Excitation and emission maxima of DsRed-Monomer

- Excitation maximum = 557 nm
- Emission maximum = 592 nm

References

1. Matz, M. V., *et al.* (1999) *Nature Biotech.* **17**(10):969–973.
2. Haas, J., *et al.* (1996) *Curr. Biol.* **6**(3):315–324.
3. Kozak, M. (1987) *Nucleic Acids Res.* **15**(20):8125–8148.
4. Gorman, C. (1985) In *DNA cloning: A Practical Approach, Vol. II*. Ed. D. M. Glover. (IRL Press, Oxford, U.K.), pp. 143–190.

Note: The attached sequence file has been compiled from information in the sequence databases, published literature, and other sources, together with partial sequences obtained by Clontech Laboratories, Inc.. This vector has not been completely sequenced.

Notice to Purchaser

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

Not-For-Profit Entities: Orders may be placed in the normal manner by contacting your local representative or Clontech Customer Service at 650.919.7300. At its discretion, Clontech grants Not-For-Profit Entities a non-exclusive, personal, limited license to use this product for non-commercial life science research use only. Such license specifically excludes the right to sell or otherwise transfer this product, its components or derivatives thereof to third parties. No modifications to the protein coding sequence may be made without express written permission from Clontech. Any other use of this product requires a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com. For-Profit Entities wishing to use this product are required to obtain a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at licensing@clontech.com or [click here](#) for more information.

The DsRed Monomer and the Fruit Fluorescent Proteins are covered by one or more of the following U.S. Patents: 7,157,566; 7,393,923; 7,005,511 and 7,250,298.

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc., unless noted otherwise. Clontech is a Takara Bio Company. ©2010 Clontech Laboratories, Inc.