


# Vector Map

## pGADT7-Rec Vector

### Catalog No.

(Not sold separately)

Sold as a part of 630490 & 630491

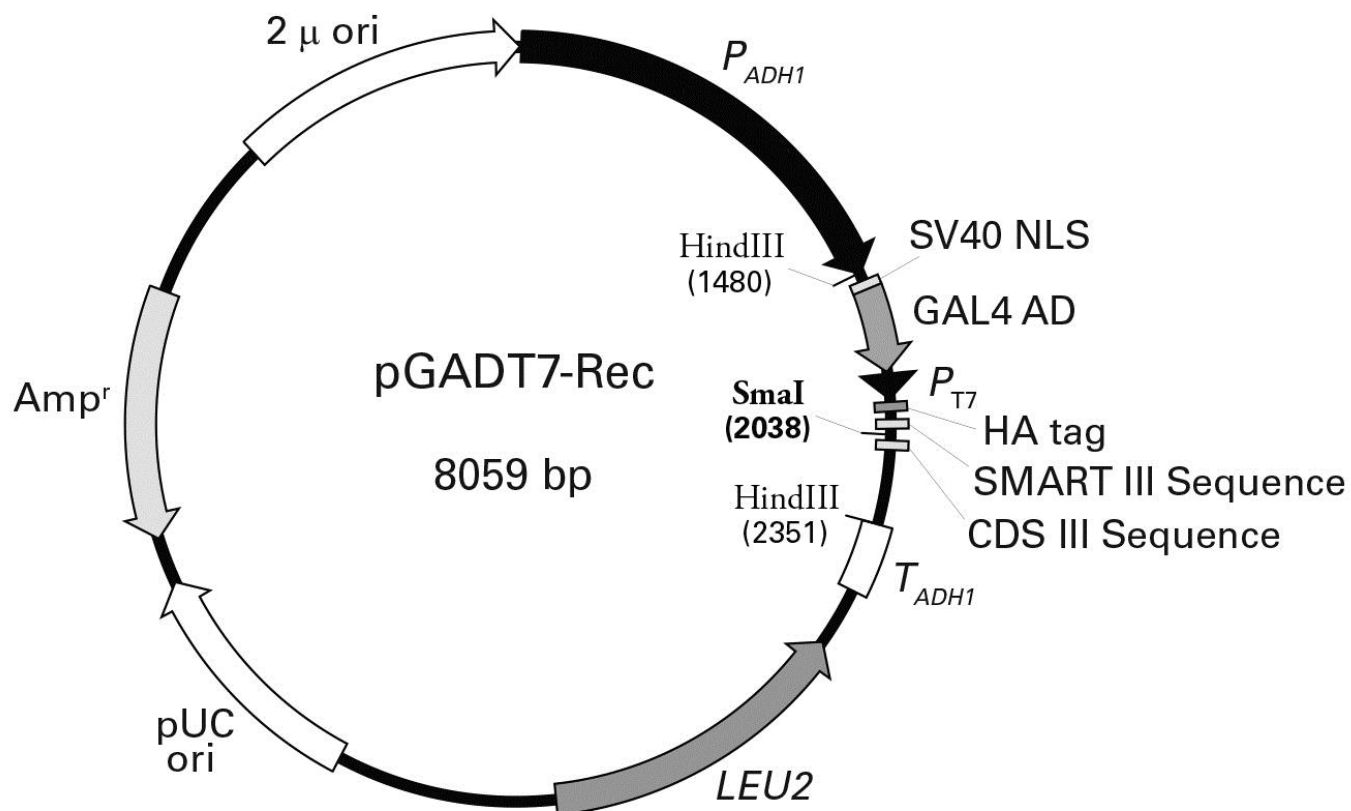

**Figure 1. pGADT7-Rec Vector map.** A unique restriction site (*Sma*I) is shown in bold. Both the Make Your Own "Mate & Plate" Library System and the Matchmaker® Gold Yeast One-Hybrid System (Cat. Nos. 630490 and 630491, respectively) contain the *Sma*I-linearized form of this vector, the form used for recombination-mediated cloning in yeast.

### Takara Bio USA, Inc.

1290 Terra Bella Avenue, Mountain View, CA 94043, USA

U.S. Technical Support: [tech@clontech.com](mailto:tech@clontech.com)


United States/Canada 800.662.2566 (102016)	Asia Pacific +1.650.919.7300	Europe +33.(0)1.3904.6880	Japan +81.(0)77.565.6999
--	---------------------------------	------------------------------	-----------------------------

# Vector Map

pGADT7-Rec Vector (Not sold separately)

Sold as a part of Cat. Nos. 630490 & 630491

## SMART III terminus

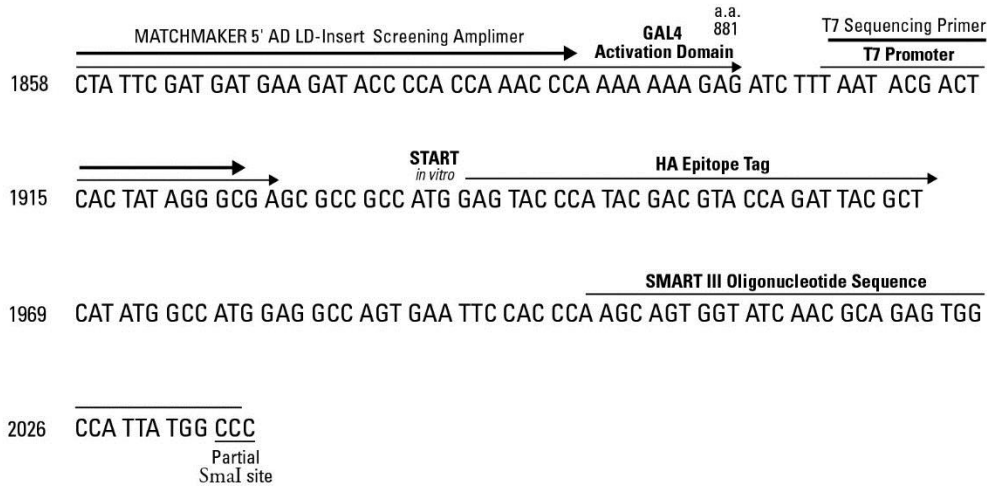

## CDS III terminus

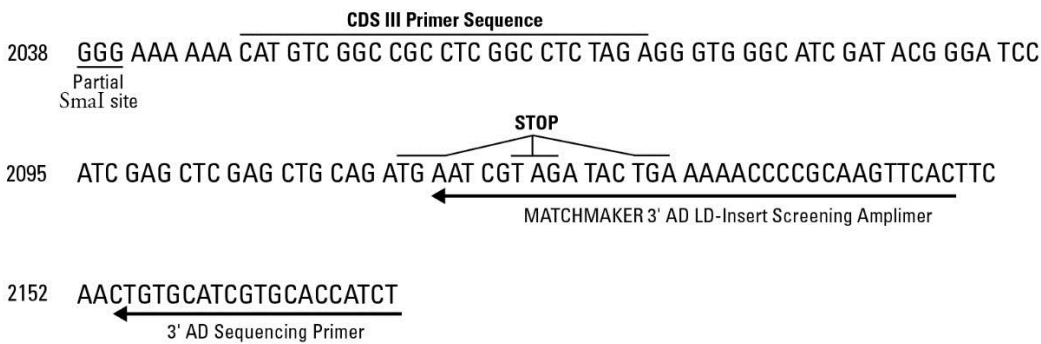


Figure 2. pGADT7-Rec Vector cloning site.

pGADT7-Rec is engineered for the construction of GAL4 AD/cDNA libraries by homologous recombination in yeast (Figure 2). Libraries made with this vector can be used for Matchmaker Gold One- and Two-Hybrid Screening.

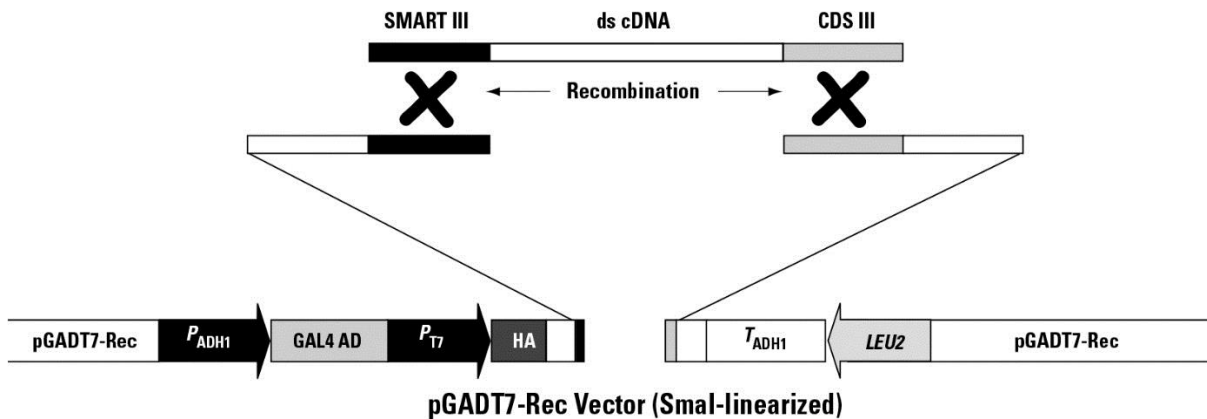


Figure 3. Cloning cDNA into pGADT7-Rec by homologous recombination *in vivo*. The ends of the SmaI-linearized vector are homologous to our SMART® III Oligonucleotide and CDS III Primer, used in the Matchmaker cDNA synthesis protocol (Figure 2).

# Vector Map

pGADT7-Rec Vector (Not sold separately)

Sold as a part of Cat. Nos. 630490 & 630491

## Location of Features

- *P<sub>ADHI</sub>* (full-length *S. cerevisiae ADHI* promoter): 7–1479
- GAL4 AD (GAL4 activation domain with SV40 nuclear localization signal [NLS]):  
SV40 NLS: 1501–1557  
GAL4 AD (amino acids 768–881): 1561–1899
- *P<sub>T7</sub>* (T7 RNA polymerase promoter): 1905–1927
- HA tag (hemagglutinin epitope tag): 1942–1968
- SMART III Oligonucleotide sequence: 2001–2036
- CDS III Primer sequence: 2047–2071
- *T<sub>ADHI</sub>* (*S. cerevisiae ADHI* Terminator): 2351–2676
- *LEU2* coding sequences: 2794–3885 (complementary)
- pUC ori (pUC replication origin): 4652–5489
- Amp<sup>R</sup> (ampicillin resistance gene): 5646–6503 (complementary)
- $\mu$  ori (yeast 2  $\mu$  replication origin): 7069–8059

## Location of Primers

- T7 Sequencing Primer: 1905–1927
- 3' AD Sequencing Primer: 2173–2154
- Matchmaker 5' AD LD-Insert Screening Amplimer (Cat. No. 630433): 1858–1889
- Matchmaker 3' AD LD-Insert Screening Amplimer (Cat. No. 630433): 2149–2117

## Propagation in *E. coli*

- Suitable host strains: DH5 $\alpha$ , DH10 & other general purpose strains
- Selectable marker: plasmid confers resistance to ampicillin (100  $\mu$ g/ml) to *E. coli* hosts
- *E. coli* replication origin: pUC
- Copy number: ~500
- Plasmid incompatibility group: pMB1/Col E1

## Propagation in *S. Cerevisiae*

- Suitable host strain: Y1HGold, Y2HGold, AH109(*MATa*), Y187(*MAT $\alpha$* ), Y190(*MATa*), SFY526(*MATa*), CG1945(*MATa*), HF7c(*MATa*)
- Selectable marker: *LEU2*
- *S. cerevisiae* origin: 2  $\mu$

**NOTE:** The vector sequence was compiled from information in the sequence databases, published literature, and other sources, together with partial sequences we obtained. This vector has not been completely sequenced.

# Vector Map

pGADT7-Rec Vector (Not sold separately)

Sold as a part of Cat. Nos. 630490 & 630491

---

## Notice to Purchaser

Our products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Our products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without prior written approval of Takara Bio USA, Inc.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product's web page at <http://www.takarabio.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

**©2016 Takara Bio Inc. All Rights Reserved.**

All trademarks are the property of Takara Bio Inc. or its affiliate(s) in the U.S. and/or other countries or their respective owners. Certain trademarks may not be registered in all jurisdictions. Additional product, intellectual property, and restricted use information is available at <http://www.takarabio.com>.

This document has been reviewed and approved by the Quality Department.